



BETA-ARRESTIN DOUBLE BRILLIANCE CONFORMATION BIOSENSORS

Summary: β -arrestins are multifunctional adaptor proteins classically associated with the “arrest” of G protein-coupled receptor (GPCR)-mediated signaling by promoting receptor desensitization and internalization. However, it is now evident that β -arrestins (mainly β -arrestin 1 and β -arrestin 2) can also scaffold signaling complexes proximal to a given GPCR and modulate the activity of various signaling networks (including ERK1/2, NF- κ B, and PI3-K). Importantly, β -arrestin-mediated signaling may be spatially and temporally distinct, and result in different biological outcomes, compared to G protein-mediated signal transduction. Consequently, GPCR-downstream signaling can occur via G protein-dependent and/or β -arrestin-dependent (G protein-independent) mechanisms (1-2). β -arrestins are recruited to activated GPCRs following receptor phosphorylation by G protein-coupled receptor kinases (GRKs) and/or other protein kinases (e.g., PKA, PKC). Different ligands acting on a given receptor are believed to produce unique receptor phosphorylation patterns (i.e., phosphorylation “barcodes”), resulting in the stabilization of different receptor-associated (active) β -arrestin conformations. Furthermore, each conformation is suggested to exhibit distinctive functional properties, thus adding an extra qualitative dimension to GPCR-downstream β -arrestin-dependent signaling (2-3).

The bioSensAll™ β -arrestin double brilliance conformation biosensors are unimolecular BRET-based biosensors that monitor conformational changes occurring within β -arrestins upon GPCR activation. These biosensors consist of β -arrestin 1 or β -arrestin 2 proteins with an N-terminal green fluorescence protein (GFP; G in following figure) tag and a C-terminal *Renilla* luciferase (RLuc; R in figure below) tag. Recruitment of β -arrestins to an activated receptor leads to a structural reorganization and subsequent physical separation of N-terminal and C-terminal regions of the β -arrestin proteins, ultimately resulting in a BRET signal decrease (4).

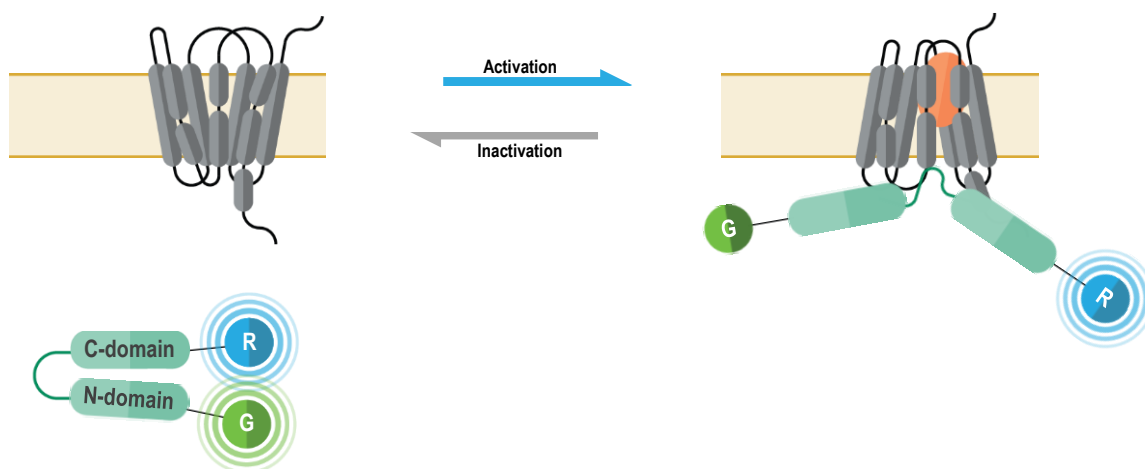

Results

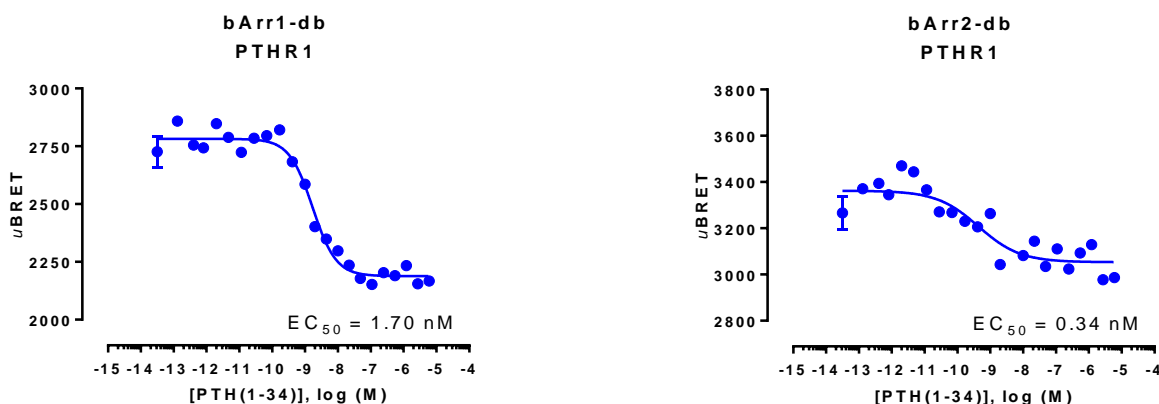

HEK293 cells were transfected with a receptor coding plasmid (human parathyroid hormone type 1 receptor (PTH1R)) in addition to the plasmid coding for the β -arrestin double brilliance conformation biosensor. On the day of BRET, cells were rinsed with assay buffer, incubated with coelenterazine and increasing amounts of PTH(1-34) for 10 minutes and BRET subsequently measured.

References

- 1- Shukla AK, Xiao K, Lefkowitz RJ. Emerging paradigms of β -arrestin-dependent seven transmembrane receptor signaling. 2011. *Trends Biochem Sci.* 36(9):457-69. PMID: 21764321.
- 2- Smith JS, Rajagopal S. The β -Arrestins: Multifunctional Regulators of G Protein-coupled Receptors. 2016. *J Biol Chem.* 291(17):8969-77. PMID: 26984408.
- 3- Yang F, Yu X, Liu C, Qu CX, Gong Z, Liu HD, Li FH, Wang HM, He DF, Yi F, Song C, Tian CL, Xiao KH, Wang JY, Sun JP. Phospho-selective mechanisms of arrestin conformations and functions revealed by unnatural amino acid incorporation and (19)F-NMR. 2015. *Nat Commun.* 6:8202. PMID: 26347956.
- 4- Zimmerman B, Beautrait A, Aguila B, Charles R, Claing A, Escher E, Bouvier M and Laporte SA. Differential β arrestin-dependent conformation signalling and cellular responses revealed by angiotensin analogs. 2012. *Sci Signal.* 5(221): ra33. PMID: 22534132.